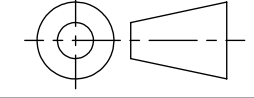



MOUNTING INSTRUCTIONS:

1. REMOVE THE FRONT BRACKET (2) FROM THE KEO SENTRY IMAGER (1) BY REMOVING THE FOUR 1/4-20 BOLTS (3).
2. MOUNT THE FRONT BRACKET (2) IN THE DESIRED LOCATION USING THE 4 1/4-20 CLEARANCE HOLES PROVIDED. THE HOLE LAYOUT OF THIS BRACKET IS SHOWN TO THE RIGHT.
3. LIFT THE KEO SENTRY IMAGER (1) UP ONTO THE MOUNTING BRACKET. IT SHOULD CLIP INTO PLACE AND BE ABLE TO HANG FREELY.
4. SECURE THE SENTRY IMAGER (1) TO THE FRONT BRACKET (2) USING THE FOUR 1/4-20 BOLTS (3) YOU PREVIOUSLY REMOVED.
5. SECURING THE REAR BRACKET (4) IS RECOMMENDED.

ITEM NO.	QTY.	DESCRIPTION
1	1	KEO ARGES-VF PHOTOMETER
2	1	FRONT MOUNTING BRACKET
3	4	1/4-20 x 1/2" SOCKET HEAD CAP SCREWS
4	1	REAR MOUNTING BRACKET

 <small>UNLESS OTHERWISE SPECIFIED DIMENSIONS ARE IN INCHES TOLERANCES ARE:</small> DECIMAL X.XX = ±0.01 X.XXX = ±0.005 DUAL DIMENSIONS 1.00in [25.4mm] REMOVE ALL BURRS AND BREAK SHARP EDGES .01 MAX.	PROPRIETARY NOTICE <small>INFORMATION CONTAINED IN THIS DOCUMENT IS THE PROPERTY OF KEO SCIENTIFIC LTD. DO NOT DISTRIBUTE OR COPY WITHOUT WRITTEN PERMISSION</small>		 Keo Scientific Ltd. <small>404 - 1300 8th St SW Calgary, AB, T2R1B2 Tel: (403) 452-7222</small>	
	APPROVALS	DATE	TITLE: Keo Arges-VF Photometer Mounting Diagram	
	DRAWN: Taylor	2012-04-23	PART NO.	REVISION
	CHECKED:		APPROVED:	C A-01
DO NOT SCALE DRAWING		SCALE = 1:2	SHEET 1 OF 1	